


Microscope Instructions

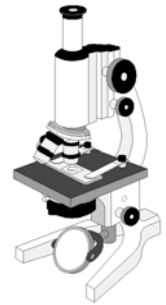

Cut out the instructions below and arrange them in order.

- a. Looking from the side, adjust the coarse focus until the objective lens is close, but not touching the slide.
- b. Select the lowest power objective lens.
- c. Use the fine focus to produce a better image.
- d. Carry the microscope carefully by the arm, with your other hand under the base.
- e. Look through the eyepiece lens and adjust the mirror so that you see a bright white circle.
- f. Look through the eyepiece and adjust the coarse focus until the image is clear.
- g. Repeat the process using a more powerful lens.
- h. Adjust the slide so that the desired part of the object is in the centre of your view.
- i. Place the microscope on the table with the arm towards you. Raise the tube as high as possible using the coarse focus.
- j. Place the slide on the stage and secure with the clips.